

FAQ

Ni Sepharose Fast Flow

HisTrap Fast Flow

HisPrep 16/10 Fast Flow



Q When should I use the new Ni Sepharose™ 6 Fast Flow instead of Ni Sepharose™ High Performance?

A The new Ni Sepharose 6 Fast Flow and Ni Sepharose High Performance are optimized for the purification of Histidine-tagged proteins and the media are pre-charged with Ni²⁺. Ni Sepharose 6 Fast Flow is designed for fast, reliable scale-up of Histidine-tagged protein purification. The medium is also well suited for batch/gravity flow purification and multi-well plate screening. Ni Sepharose High Performance is optimized for high performance purification of Histidine-tagged proteins with a chromatography system, such as ÄKTAdesign from Amersham Biosciences.

Q What is the protein binding capacity of Ni Sepharose 6 Fast Flow?

A *Dynamic binding capacity:* Approximately 40 mg (His)₆-tagged protein/ml medium

Dynamic binding capacity conditions:

<i>Sample:</i>	1 mg/ml (His) ₆ -tagged pure protein (M _r 43 000) in binding buffer (QB _{10%} determination) or (His) ₆ -tagged protein (M _r 28 000) bound from <i>E. coli</i> extract
<i>Column volume:</i>	0.25 ml or 1 ml
<i>Flow rate:</i>	0.25 ml/min or 1 ml/min
<i>Binding buffer:</i>	20 mM sodium phosphate, 0.5 M NaCl, 5 mM imidazole, pH 7.4
<i>Elution buffer:</i>	20 mM sodium phosphate, 0.5 M NaCl, 500 mM imidazole, pH 7.4

Note: Dynamic binding capacity is protein dependent.

Q What are the general running conditions recommended?

A **Recommended starting conditions:**

<i>Binding buffer:</i>	20 mM sodium phosphate, 0.5 M NaCl, 20–40 mM imidazole, pH 7.4
<i>Elution buffer:</i>	20 mM sodium phosphate, 0.5 M NaCl, 500 mM imidazole, pH 7.4

Note: The concentration of imidazole that will give optimal purification results (in terms of purity and yield) is protein-dependent, and is usually slightly higher for Ni Sepharose 6 Fast Flow than for similar IMAC media on the market. This can be useful during the washing step as a higher concentration of imidazole in the wash buffer may wash out contaminants that have bound to the medium. This can result in elution of a more pure target protein.

Higher concentration of imidazole will not cause any problems in terms of increased baseline absorbance if a high purity imidazole is used, such imidazole gives essentially no absorbance at 280 nm.

Q Will the high capacity of the new Ni Sepharose 6 Fast Flow result in higher non-specific binding?

A The non-specific binding of proteins is a result of several factors. Other proteins will only bind more to Ni Sepharose 6 Fast Flow if these non-specific proteins are interacting directly with the Ni²⁺ or the metal chelating ligand. In addition, non-specific proteins can potentially be washed off at higher concentrations of imidazole, while keeping the target protein bound to the medium.

Q With the higher binding capacity, do you have to use harsh conditions to remove the target protein?

A Harsh conditions are not necessary to remove most target proteins. The proteins tested are all eluted below 500 mM imidazole; frequently between 100-300 mM imidazole. The protein is efficiently eluted from the column, resulting in high yields.

Q How many times can you run the Ni Sepharose 6 Fast Flow while maintaining the same capacity?

A This is sample dependent. Reusing Ni Sepharose 6 Fast Flow depends on the nature of the sample. To prevent cross-contamination, only reuse the column when purifying identical Histidine-tagged proteins.

Q How do you measure the amount of Ni²⁺ in the eluent?

A This quantitative measurement assay is as follows:

The medium is run with 10 column volumes of a buffer with pH 4.0 and then the Ni²⁺ still bound on the column is measured spectroscopically and compared with the amount Ni²⁺ loaded.

Q Can you strip and regenerate the Ni Sepharose 6 Fast Flow medium?

A The new Ni Sepharose 6 Fast Flow can be stripped with 50 mM EDTA in 20 mM sodium phosphate, 0.5 M NaCl, pH 7.4 and recharged with Ni²⁺ or other divalent cations, such as Cu²⁺, Co²⁺, Zn²⁺, etc.

Q How can I strip and recharge the medium?

A Recommended stripping buffer: 20 mM sodium phosphate, 0.5 M NaCl, 50 mM EDTA, pH 7.4.

1. Strip the column by washing with 5–10 column volumes of stripping buffer.
2. Wash with 5–10 column volumes of binding buffer.
3. Wash with 5–10 column volumes of distilled water.
4. Re-charge the column.

To recharge the water-washed column:

1. Load 0.5 column volumes of 0.1 M metal salt solution in distilled water on the column. Metal chloride and sulfate salts, e.g. 0.1 M NiSO₄, are commonly used.
2. Wash with 5 column volumes of distilled water.
3. Wash with 5 column volumes of binding buffer.

Note: It is important to wash with binding buffer as the last step to get pH correct.

Q How is the Ni²⁺ leakage at low pH?

A At pH 4, the loss of Ni²⁺ is < 5% on Ni Sepharose 6 Fast Flow (reproducible) compared to average 9% for Ni-NTA Superflow from Qiagen (wide batch to batch variations).

Note: Loss of metal ions is more pronounced at lower pH.

Q My target protein is extremely sensitive to Ni²⁺. How do I reduce the Ni²⁺ leakage to a real minimum (below <5 %)?

A Leakage of Ni²⁺ from Ni Sepharose 6 fast Flow is low under all conditions and resistance towards reducing agents is thus high.

Note: If using buffers/sample including reducing agents it is recommended to perform a blank run *without* reducing agents in the buffers before loading sample (to remove any weakly bound Ni²⁺ ions):

Blank run: Use binding and elution buffer *without* reducing agents

1. Wash the column with 5 column volumes of distilled water.
2. Wash with 5 column volumes of binding buffer.
3. Wash with 5 column volumes of elution buffer.
4. Equilibrate with 10 column volumes of binding buffer.

Q What different additives can I use together with Ni Sepharose 6 Fast Flow, HisPrep™ FF 16/10 and HisTrap™ FF?

A Tests performed in our laboratories show that Ni Sepharose 6 fast Flow is compatible with the following compounds at the concentrations given.

Reducing agents	5 mM DTE	
	5 mM DTT	
	20 mM β-mercaptoethanol	
	5 mM TCEP	
	10 mM reduced glutathione	
Denaturing agents	8 M urea*	
	6 M guanidine hydrochloride*	
Detergents	2% Triton™ X-100	Non-ionic detergent
	2% Tween™ 20	Non-ionic detergent
	2% NP-40	Non-ionic detergent
	2% cholate	Anionic Detergent
	1% CHAPS	Zwitterionic Detergent
Additives	500 mM imidazole	
	20% ethanol	
	50% glycerol	
	100 mM Na ₂ SO ₄	
	1.5 M NaCl	
	1 mM EDTA**	
	60 mM citrate**	
Buffer solutions	50 mM sodium phosphate, pH 7.4	
	100 mM Tris-HCl, pH 7.4	
	100 mM Tris-acetate, pH 7.4	
	100 mM HEPES, pH 7.4	
	100 mM MOPS, pH 7.4	
	100 mM sodium acetate, pH 4*	

* Tested for 1 week at +40 °C

** Generally, chelating agent should be used with caution (and only in the sample, not in buffers). Any metal ion stripping may be counteracted by addition of a small excess of MgCl₂ before centrifugation or filtration of the sample. Note that stripping effects may vary with applied sample volume.

Q How do I remove imidazole from my target protein after elution?

A If imidazole needs to be removed from the protein, use HiTrap™ Desalting 5 ml (can be coupled in series to increase the sample volume), a PD-10 column, or HiPrep™ 26/10 Desalting, depending on the sample volume.

Q Do you see an improved purity (less non-specific binding) with high concentrations of NaCl?

A In most cases, concentration of ions, imidazole, and other additives in the binding and wash buffers will influence the level of non-specific binding. Higher concentrations of NaCl (500 mM) generally reduce the level of non-specific binding.

Q How does DTT affect the performance of Ni Sepharose 6 Fast Flow in comparison to nickel media from other tested suppliers?

A DTT has a drastic effect on the colour and performance of nickel medium from other tested suppliers, so DTT is generally avoided.

For Ni Sepharose 6 Fast Flow, however, DTT has been tested up to 5 mM and does not affect the binding capacity or nickel ion leakage of the Ni Sepharose 6 Fast Flow. There is no effect on the performance of the medium, although color changes can be evident over 3mM DTT. *Therefore, it is essential to perform a blank run before loading sample/buffers including DTT (and other reducing agents).*

Blank run: Use binding and elution buffer *without* reducing agents

1. Wash the column with 5 column volumes of distilled water.
2. Wash with 5 column volumes of binding buffer.
3. Wash with 5 column volumes of elution buffer.
4. Equilibrate with 10 column volumes of binding buffer.

Q Is it possible to use high concentrations of Tris-HCl buffer and how does this compatibility to Tris buffers compare to other suppliers?

A These are the Tris buffers systems that have been tested and work well.

50 mM Tris-HCl, pH 7.4

100 mM Tris-HCl, pH 7.4

100 mM Tris-acetate, pH 7.4

Tris buffers in high concentrations can adversely affect the binding of Histidine-tagged target proteins to nickel medium from other suppliers; therefore, lower concentrations of Tris buffers (less than 50 mM) or phosphate buffers are generally used.

Q How can I clean Ni Sepharose 6 Fast Flow?

A Note: Strip off the metal ions before cleaning (to prevent precipitation of metal salt)

- To remove precipitated proteins, hydrophobically bound proteins and lipoproteins:
 1. Wash the column with 1 M NaOH, contact time usually 1–2 hours (12 hours or more for endotoxin removal).
 2. Wash the column with approximately 10 column volumes of binding buffer.
 3. Wash the column with 5–10 column volumes of distilled water.
- To remove strongly hydrophobically bound proteins, lipoproteins and lipids:
 1. Wash the column with 5–10 column volumes 30% isopropanol for about 15–20 minutes.
 2. Wash the column with approximately 10 column volumes of distilled water.
 3. Re-charge the medium or wash with 5 column volumes 20% ethanol for storage.

Alternatively:

1. Wash the column with 2 column volumes of detergent in a basic or acidic solution. Use, for example, 0.1–0.5% nonionic detergent in 0.1 M acetic acid, contact time 1–2 hours.
 2. After treatment, always remove residual detergent by washing with 5 column volumes of 70% ethanol.
 3. Wash the column with approximately 10 column volumes of distilled water.
 4. Re-charge the medium or wash with 5 column volumes 20% ethanol for storage.
- To remove ionically bound proteins:
 1. Washing the column with 1.5 M NaCl solution, contact time 10–15 minutes.
 2. Wash the column with approximately 10 column volumes of distilled water.
 3. Re-charge the medium or wash with 5 column volumes 20% ethanol for storage.

Q Is it possible to use the new Ni Sepharose 6 Fast Flow by gravity or in 96 or 384 well multi-well plates for screening?

A Yes, the performance by gravity flow has been tested. The new Ni Sepharose 6 Fast Flow can be used to fill empty gravity columns and multi-well plates.

Q Do the new HisTrap FF columns have Valco fittings?

A Yes. This makes it possible to put these columns directly into ÄKTAdesign™ without any extra connectors.

www.gehealthcare.com

GE Healthcare
Amersham Biosciences AB
Björkgatan 30
751 84 Uppsala
Sweden

ÅKTAdesign, HisPrep, HisTrap and Sepharose are trademarks of GE Healthcare Ltd, a General Electric Company. GE Tagline and GE Monogram are trademarks of General Electric Company.

Triton is a trademark of Union Carbide Chemicals and Plastics Co. Tween is a trademark of ICI Americas Inc. Qiagen is a trademark of the Qiagen group. Superflow is a trademark of Sterogene Bioseparations Inc.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. GE Healthcare reserves the right, subject to any regulatory and contractual approval, if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation.

© 2005 General Electric Company – All rights reserved.

Amersham Biosciences AB, a General Electric Company going to market as GE Healthcare.

GE Healthcare Amersham Biosciences AB
Björkgatan 30, 751 84 Uppsala, Sweden

GE Healthcare Amersham Biosciences Europe GmbH
Munzinger Strasse 9, D-79111 Freiburg, Germany

GE Healthcare Amersham Biosciences UK Ltd
Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare Amersham Biosciences Corp
800 Centennial Avenue, P.O. Box 1327, Piscataway, NJ 08855-1327, USA

GE Healthcare Amersham Biosciences KK
Sanken Bldg. 3-25-1, Hyakunincho, Shinjuku-ku, Tokyo 169-0073, Japan

GE Healthcare Amersham Biosciences Ltd
13/F, Tower 1, Ever Gain Plaza, 88 Container Port Road Kwai Chung, New Territories, HONG KONG

Asia Pacific Tel: +852 2811 8693 Fax: +852 2811 5251 • Australasia Tel: + 61 2 9899 0999 Fax: +61 2 9899 7511 • Austria Tel: 01/57606-1619 Fax: 01/57606-1627 • Belgium Tel: 0800 73 888 Fax: 03 272 1637 • Canada Tel: 800 463 5800 Fax: 800 567 1008 • Central, East, & South East Europe Tel: +43 1 982 3826 Fax: +43 1 985 8327 • Denmark Tel: 45 16 2400 Fax: 45 16 2424 • Finland & Baltics Tel: +358-09-512 39 40 Fax: +358 (0)9 512 39 439 • France Tel: 01 69 35 67 00 Fax: 01 69 41 96 77 • Germany Tel: 0761/4903-490 Fax: 0761/4903-405 • Italy Tel: 02 27322 1 Fax: 02 27302 212 • Japan Tel: +81 3 5331 9336 Fax: +81 3 5331 9370 • Latin America Tel: +55 11 3933 7300 Fax: +55 11 3933 7304 • Middle East & Africa Tel: +30 210 9600 687 Fax: +30 210 9600 693 • Netherlands Tel: 0165 580 410 Fax: 0165 580 401 • Norway Tel: 815 65 555 Fax: 815 65 666 • Portugal Tel: 21 417 7035 Fax: 21 417 3184 • Russia & other C.I.S. & N.I.S Tel: +7 (095) 232 0250 Fax: +7 (095) 230 6377 • South East Asia Tel: 60 3 8024 2080 Fax: 60 3 8024 2090 • Spain Tel: 93 594 49 50 Fax: 93 594 49 55 • Sweden Tel: 018 612 1900 Fax: 018 612 1910 • Switzerland Tel: 0848 8028 12 Fax: 0848 8028 13 • UK Tel: 0800 616928 Fax: 0800 616927 • USA Tel: 800 526 3593 Fax: 877 295 8102

