

# illustra AutoSeq G-50 and AutoSeq 96 G-50

## Introduction

illustra™ AutoSeq™ G-50 and the new illustra AutoSeq 96 G-50 products provide a convenient and simple solution for dye terminator removal from sequencing samples prior to analysis. illustra AutoSeq G-50 is useful for handling small numbers of samples and illustra AutoSeq 96 G-50 is designed for parallel processing of up to 96 samples (Fig 1). We introduce illustra AutoSeq 96 G-50 for use with MegaBACE™ and illustra AutoSeq 96 G-50 for use with Applied Biosystems Inc., (ABI) sequence analyzers, respectively.

Purification of large numbers of sequencing reactions using single columns is time consuming, expensive, and tedious. In addition, ethanol precipitation—which is the standard laboratory method for dye terminator removal—is inconsistent with regards to sample recovery and purity. illustra AutoSeq G-50 and AutoSeq 96 G-50 provide faster, more consistent results than ethanol precipitation, with greater ease-of-use for multiple samples.

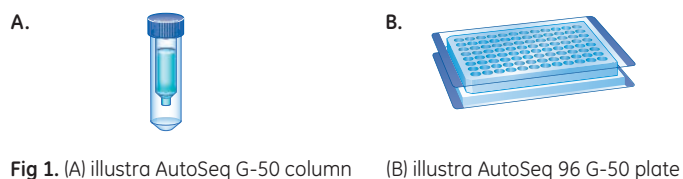


Fig 1. (A) illustra AutoSeq G-50 column (B) illustra AutoSeq 96 G-50 plate

### illustra AutoSeq G-50 and AutoSeq 96 G-50 deliver:

- **Speed:** Pre-equilibrated columns and 96-well plates allow you to go from sample application to collection of purified DNA product in less than 4 min for the column and 15 min for the 96-well plate
- **Simplicity:** Quick reference protocol card provides instructions at a glance for experienced users
- **High-quality DNA:** Provides equivalent Phred 20 read length results as compared with DyeEx™ 2.0 Spin Kit from Qiagen™ and Edge Biosystems' Performa™ V3 DTR 96-Well Short Plate

## Method overview

The use of relatively small volumes (5 or 10  $\mu$ l) in sequencing reactions is becoming increasingly popular with researchers who perform high-throughput sequencing on ABI analyzers because of reduced costs while maintaining good signal strength and Phred 20 read length. illustra AutoSeq products contain a pre-equilibrated and pre-weighed amount of Sephadex™ G-50 DNA media for optimal sequencing reaction purification. The purification process involves gel filtration and centrifugation to separate sequencing reaction products from excess dye terminators. The exclusion properties that allow the sequencing product to pass through the gel whilst retaining the smaller dye terminator and salt impurities depend on factors such as the type of resin used, sample volume, product size, and the  $g$  forces employed in the purification process. The protocols provided with the illustra AutoSeq products have been optimized for the purification of dye terminator sequencing reactions (Fig 2).

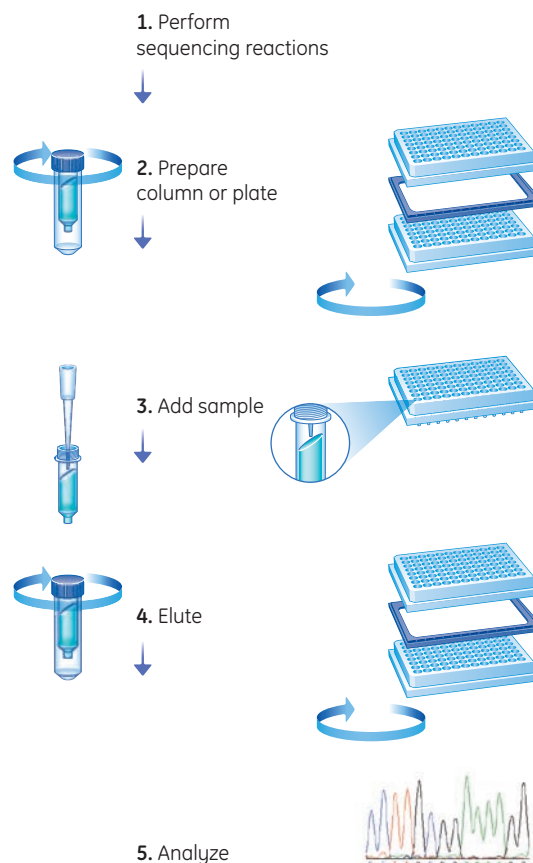



Fig 2. Schematic representation of the method employed by illustra AutoSeq G-50 dye terminator removal products.



## High-quality sequencing analysis

Phred quality (*q*) values express the probability of correctly calling a base-peak in a sequencing electropherogram and they are based on key parameters such as resolution, spectral cross talk, and uniformity of peak spacing (1–3). The *q* value is logarithmically proportional to the probability of correctly calling the identified peak (i.e., a Phred value of 20 is an accuracy of 99%, Phred 30 is 99.9% accurate). The minimal acceptable value for sequencing is a Phred value of 20. For this study, we defined sequencing read length as the total number of bases with a Phred value of 20 or greater ( $q > 20$ ) using a 10-base moving window average.

Tests using 20  $\mu$ l reaction volumes showed no significant differences in Phred 20 read length obtained with the Qiagen DyeEx 2.0 Spin Kit and the illustra AutoSeq G-50 dye terminator removal kit (Fig 3) indicating that the illustra AutoSeq G-50 dye terminator removal kit can be used to generate high-quality sequencing results.

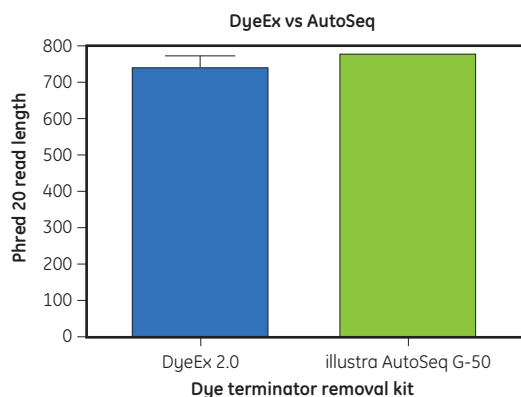


**Fig 3.** Phred 20 read length obtained using DyeEx 2.0 and AutoSeq G-50 kits ( $p = 0.2061$ ). A bulk stock of sequencing reactions with BigDye™ Terminator v3.1 Cycle Sequencing Kit was prepared using a pUC-based plasmid containing the TP53 open reading frame (Tumor protein p53) as template and M13-forward or -reverse sequencing primers. Aliquots from this bulk stock were purified. Three sequencing reactions using the forward primer and three reactions using the reverse primer were purified by 3 different researchers. Purified sequencing samples were analyzed by the Qiagen Sequencing Service for analysis using an ABI PRISM™ 3730XL DNA analyzer.


We checked for data accuracy at the beginning of the sequence by examining electropherograms for the absence of dye blobs and noting the read length at which the Phred score reached  $\geq 20$ . Dye blobs were absent from all the samples tested and there was no significant difference in the read length at which the Phred score reached 20 or more ( $p = 0.6579$ ) between illustra AutoSeq-G50 and DyeEx 2.0 columns. A typical electropherogram image depicting the quality and accuracy of data generated with the illustra AutoSeq-G50 Kit is shown in (Fig 4).

## High-throughput sequencing analysis

Test reactions using both 5 and 10  $\mu$ l sequencing reaction volumes and 1  $\mu$ l of dye terminator mix were performed. There was no significant difference in the Phred 20 read lengths for reactions purified with the illustra AutoSeq 96 G-50, Qiagen's DyeEx 96 Kit, and Edge Biosystems' Performa product (Fig 5).

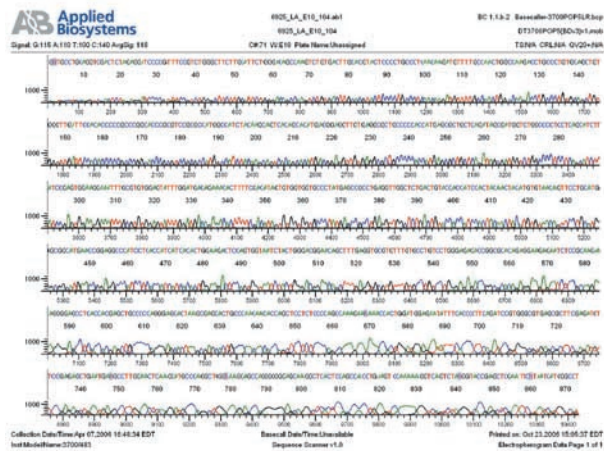

**Fig 4.** Typical electropherogram of sequencing reactions purified using the illustra AutoSeq G-50.

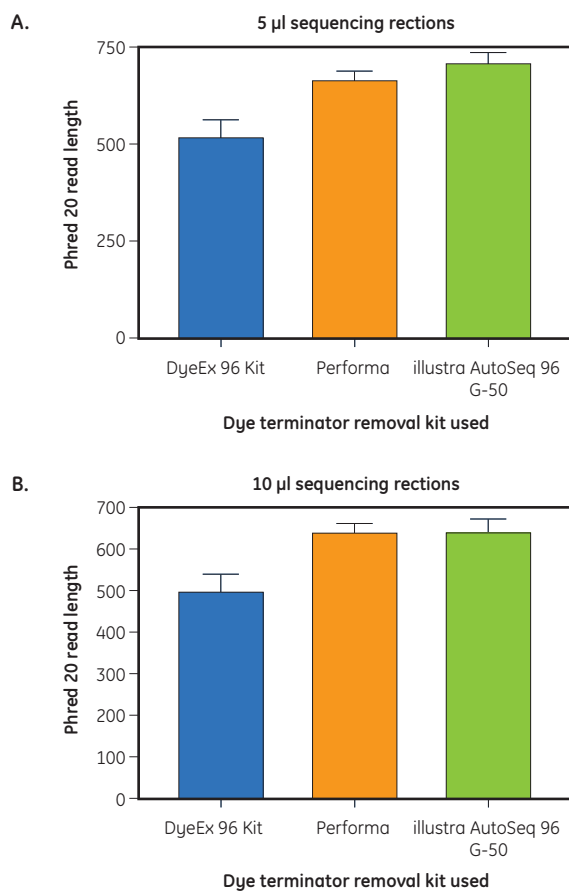

**Fig 5.** Phred 20 read length obtained from different reaction volumes ( $p > 0.05$ ). A bulk stock of sequencing reactions was prepared with ABI's BigDye Terminator v3.1 Cycle Sequencing Kit using a pUC-based plasmid containing the Tumor protein p53 open reading frame (ORF) as template and M13-forward or -reverse sequencing primers. Aliquots from this bulk stock were purified. Purified sequencing samples were analyzed by the Qiagen Sequencing Service for analysis on an ABI 3730XL. To test for well-to-well variability, 6 samples of forward primers and another 6 samples of reverse primers were purified with the 96-well purification kits.

## Summary

illustra AutoSeq G-50 and illustra AutoSeq 96 G-50 kits are designed for the rapid removal of unincorporated dye terminators from sequencing reactions. The purified DNA product routinely generates high-quality sequencing results. We now provide two new AutoSeq 96-well format kits for use on MegaBACE and ABI sequence analyzers, respectively. The versatility of the illustra AutoSeq product range enables you to take advantage of using small volumes of reagents for sequencing reactions without compromising sequence quality.

### References:

1. Ewing, B. and Green, P. Base-calling of automated sequencer traces using phred. II. Error probabilities. *Genome Res.* **8**, 186-194 (1998).
2. Ewing, B *et al.* Base-calling of automated sequencer traces using phred. I. Accuracy assessment. *Genome Res.* **8**, 175-185 (1998).
3. Paegel, B. M. *et al.* High throughput DNA sequencing with a microfabricated 96-lane capillary array electrophoresis bioprocessor. *Proc. Natl. Acad. Sci. U S A.* **99**, 574-579. (2002).

## Ordering information

illustra AutoSeq G-50 (50 columns)	27-5340-01
illustra AutoSeq G-50 (250 columns)	27-5340-02
illustra AutoSeq G-50 (1000 columns)	27-5340-03
illustra AutoSeq 96 G-50 for MegaBACE Analyzers (10 × 96-well plates)	28-9094-90
illustra AutoSeq 96 G-50 for ABI Analyzers (10 × 96-well plates)	28-9034-27

**Asia Pacific** Tel: +65 6275 1830 Fax: +65 6275 1829 **Australasia** Tel: + 61 2 9899 0999 Fax: +61 2 9899 7511 **Austria** Tel: 01/57606-1619 Fax: 01/57606-1627  
**Belgium** Tel: 0800 73 888 Fax: 02 416 82 06 **Canada** Tel: 1 800 463 5800 Fax: 1 800 567 1008 **Central, East, & South East Europe** Tel +43 1 972720  
Fax: +43 1 97272 2750 **Denmark** Tel: 45 16 2400 Fax: 45 16 2424 **Eire** Tel: 01494 544000 Fax: 0044 1494 542010 **Finland & Baltics** Tel: +358-(0)9-512 39 40  
Fax: +358 (0)9 512 39 439 **France** Tel: 01 6935 6700 Fax: 01 6941 9677 **Germany** Tel: (089) 96281 660 Fax: (089) 96281 620 **Greater China** Tel: +852 2100 6300  
Fax: +852 2100 6338 **Italy** Tel: 02 27322 1 Fax: 02 27302 212 **Japan** Tel: +81 3 5331 9336 Fax: +81 3 5331 9370 **Latin America** Tel: +55 11 3933 7300  
Fax: +55 11 3933 7304 **Middle East & Africa** Tel: +30 210 9600 687 Fax: +30 210 9600 693 **Netherlands** Tel: 0800 82 82 82 1 Fax: 0800 82 82 82 4  
**Norway** Tel: 815 65 555 Fax: 815 65 666 **Portugal** Tel: 21 417 7035 Fax: 21 417 3184 **Russia & other C.I.S. & N.I.S** Tel: +7 (495) 956 5177 Fax: +7 (495) 956 5176  
**Spain** Tel: 93 594 49 50 Fax: 93 594 49 55 **Sweden** Tel: 018 612 1990 Fax: 018 612 1910 **Switzerland** Tel: 0848 8028 12 Fax: 0848 8028 13 **UK** Tel: 0800 616928  
Fax: 0800 616927 **USA** Tel: +1 800 526 3593 Fax: +1 877 295 8102

[www.gelifesciences.com/illustra](http://www.gelifesciences.com/illustra)

GE Healthcare UK Limited  
Amersham Place  
Little Chalfont  
Buckinghamshire  
HP7 9NA  
UK



imagination at work

GE, imagination at work and GE monogram are trademarks of General Electric Company.

AutoSeq, illustra, Sephadex, and MegaBACE are trademarks of GE Healthcare companies.

DyeEx and Qiagen are trademarks of the Qiagen Group. ABI PRISM and BigDye are trademarks of Applied Biosystems Inc. Performa is a trademark of Edge Biosystems.

© 2007 General Electric Company - All rights reserved. First published March 2007

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request.

Contact your local GE Healthcare representative for the most current information.

GE Healthcare Bio-Sciences AB  
Björkgatan 30, 751 84 Uppsala  
Sweden

GE Healthcare Europe, GmbH  
Munzinger Strasse 5, D-79111 Freiburg  
Germany

GE Healthcare Bio-Sciences Corp.  
800 Centennial Avenue, P.O. Box 1327  
Piscataway, NJ 08855-1327  
USA

GE Healthcare Bio-Sciences KK  
Sanken Bldg., 3-25-1, Hyakunincho  
Shinjuku-ku, Tokyo 169-0073  
Japan