


Series S Sensor Chip Protein G

A Biacore Extend product

This product is to be considered a development product and not a standard Biacore consumables product. For more information, see www.gelifesciences.com/biacoreextend.

Product description

Order code:	29179315 (package of one sensor chip)
Storage:	The use-before date applies to chips stored at 2°C to 8°C in unopened pouches.

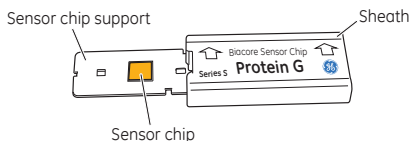

The sensor chip is fixed to a polystyrene support sheath. Each cassette, consisting of a sensor chip and sheath assembly, is individually packed under a nitrogen atmosphere in a sealed pouch.

Sensor Chip Protein G consists of a carboxymethylated dextran matrix pre-immobilized with a recombinant Protein G - GammaBind G, Type 2.

Note: For *in vitro* use only.



Application areas

Series S Sensor Chip Protein G is designed to bind antibodies of different species and subclasses for interaction analysis in Biacore systems. Series S Sensor Chip Protein G is a good choice for kinetic characterization and concentration analyses in a wide range of applications.

Refer to www.gelifesciences.com/biacore for updates on applications and scientific publications, and to *Biacore Assay Handbook* (29019400).

Surface specificity

The recombinant Protein G binds a broad range of IgG, such as human (including IgG₃), rat, rabbit, mouse, guinea pig, goat, sheep and cow.

Preparations for use

Step	Action
1	Allow the sealed sensor chip pouch to equilibrate at room temperature for 15 to 30 minutes in order to prevent condensation on the chip surface.
2	Prepare the Biacore instrument with running buffer. The buffer should be filtered (0.22 µm), and degassed for systems that do not have an integrated buffer degasser.
3	Open the sensor chip pouch. Make sure that the sensor chip support remains fully inserted into the sheath at all times to protect the chip from dust particles.
4	Dock the sensor chip in the instrument as described in the instrument handbook.

Note: *Storage stability is affected by exposure to air. Keep sensor chip in unopened pouch until use.*

Analysis temperature

Series S Sensor Chip Protein G is designed for use at 25°C.

Start-up cycles

For best assay performance, run at least one start-up cycle using sample or buffer as analyte and identical settings as for the analysis cycles.

Regeneration

Regenerate the surface with one 30-second injection of 10 mM Glycine-HCl, pH 1.5 (available from GE, product code BR100354). This will remove captured ligand together with any analyte bound to them.

Alternative regeneration procedures are:

- One additional 30 s injection of 0,05-0,5% Sodium dodecyl sulfate (SDS), or
- One 30 s injection of 10 mM Glycine-HCl pH 1.5 with 0,5% Surfactant P20 added (available from GE, product code BR100054).

Avoid using basic regeneration solutions with a pH >10

For local office contact information, visit
www.gelifesciences.com/contact

GE Healthcare UK Limited
Amersham Place
Little Chalfont
Buckinghamshire, HP7 9NA
United Kingdom

www.gelifesciences.com/biacore

GE, GE monogram and Biacore are trademarks of General Electric Company.

© 2016 General Electric Company.
First published Jan. 2016.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

GE Healthcare Bio-Sciences AB
Björkgatan 30, 751 84 Uppsala, Sweden

GE Healthcare Europe GmbH
Munzinger Strasse 5, D-79111 Freiburg, Germany

GE Healthcare Bio-Sciences Corp.
100 Results Way, Marlborough, MA 01752, USA

GE Healthcare Dharmacon, Inc.
2650 Crescent Dr., Lafayette, CO 80026, USA

HyClone Laboratories, Inc.
925 W 1800 S, Logan, UT 84321, USA

GE Healthcare Japan Corporation
Sanken Bldg. 3-25-1, Hyakunincho Shinjuku-ku, Tokyo 169-0073, Japan

